
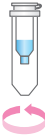

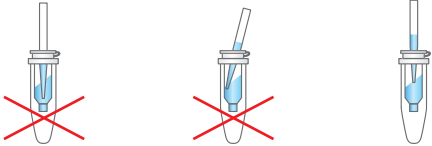



FastGene™ Dye Terminator Removal kit

For Removal of Sequencing Dye Terminator

FG-9411 50preps

Resin preparation	 <p>6.3ml Buffer DT Vortexing >30min at room temperature</p>
Column preparation	 <p>Apply 600 µl hydrated resin into the column 750 x g ; 3min</p>
	 <p>Transfer the column into a new tube.</p>
Loading sample	 <p>Up to 20 µl of sample solution</p> <p>Do not touch the gel surface with the tip.</p> <p>Do not load the sample on the inner wall of the column.</p> <p>Load the sample gently onto the center of the gel.</p>
Sample purification	 <p>750 x g ; 3min</p>

10P1105DT