

FAS-V

Quick Guide v 1.0



Lens adjustments

APERTURE RING

- Start at 4 for best sharpness/brightness
 - 1.2 for brightening the picture (for dark gels)
 - 16 for darker pictures (sharper images)

ZOOM RING

- Start at 12.5 mm (all imaging area is covered)
- Zoom to 75 mm for single bands

FOCUSSING RING

- Set focus once
- There is no need to adjust it afterwards



Taking an image

1. Set the aperture ring to 4
2. Adjust the exposure time by pressing the **EXPOSURE** button
 - a. Start at 1.5 seconds
 - b. Reduce if image is too bright
 - c. Increase if image is too dark
3. Press the CAPTURE button to take an image.
4. Choose one of the 6 precaptured images
5. Save the image