



## Restriction Enzyme Msp I



Cat.#	Size	Conc.
FG-Mspl	5,000 units	20 units/µl

Store at -20°C

**Supplied with:** 10X FastGene® Buffer IV (FG-REB4)  
10X FastGene® FastCut Buffer (FG-REBHF)  
6X DNA Loading Buffer  
Sterile water

### Recognition site



For Research Use Only. Not for use in diagnostic procedures.

ISO9001

**Source:** *Moraxella* species

### Reaction conditions

1X FastGene® Buffer IV 37°C  
1X FastGene® FastCut Buffer, 37°C

### FastGene® FastCut Buffer

FastGene® restriction enzyme can cut substrate DNA in 5-15 with FastGene® FastCut Buffer.

### 1X FastGene® Buffer IV

20 mM Tris-acetate (pH 7.9 at 25°C)  
50 mM potassium acetate  
10 mM magnesium acetate  
100 µg/ml BSA

### Unit definition

One unit is defined as the amount of enzyme required for complete digestion of 1 µg bacteriophage λ at 37°C for 1 hr in 50 µl reaction mixtures.

### Quality control

- Unit definition assay
- Overdigestion assay
- Endonuclease assay
- Extreme pure assay

### Dilution buffer:

FastGene® Diluent A

### Heat Inactivation

No.

### Methylation sensitivity

*dam* methylation: Not sensitive  
*dcm* methylation: Not sensitive  
CpG methylation: Not sensitive

### Prolonged incubation

A minimum amount of enzyme required to digest 1 µg substrate DNA for 16 hr; 0.5 U.

### Relative activity in FastGene® Buffers

FastGene® Buffer I: 75%  
FastGene® Buffer II: 100%  
FastGene® Buffer III: 75%  
FastGene® Buffer IV: 100%  
FastGene® FastCut Buffer: 100%

### Note

It is an isoschizomer of Hpa II. It is not affected by *dam*, *dcm*, or mammalian CpG methylation. However MspI can not digestion, when the external C in the sequence CCGG is methylated.

### Standard reaction condition

- Normal protocol

Component	Final Conc.	Volume
Substrate DNA	1 µg	X µl
10X FastGene® Buffer IV	1 X	5 µl
Msp I	20 unit	1 µl
Sterile water		up to 50 µl

→ Incubate at 37°C for 1 hr

- Fast protocol

Component	Final Conc.	Volume
Substrate DNA	1 µg	X µl
10X FastGene® FastCut Buffer	1 X	5 µl
Msp I	20 unit	1 µl
Sterile water		up to 50 µl

→ Incubate at 37°C for 15 min

※'Standard conditions' is only a general recommendation. The experimental conditions should be adjusted according to the purpose and sample.



## Restriction Enzyme Msp I



Cat.#	Size	Conc.
FG-Mspl	5,000 units	20 units/µl

Store at -20°C

**Supplied with:** 10X FastGene® Buffer IV (FG-REB4)  
10X FastGene® FastCut Buffer (FG-REBHF)  
6X DNA Loading Buffer  
Sterile water

### Recognition site



For Research Use Only. Not for use in diagnostic procedures.

ISO9001

**Source:** *Moraxella* species

### Reaction conditions

1X FastGene® Buffer IV 37°C  
1X FastGene® FastCut Buffer, 37°C

### FastGene® FastCut Buffer

FastGene® restriction enzyme can cut substrate DNA in 5-15 with FastGene® FastCut Buffer.

### 1X FastGene® Buffer IV

20 mM Tris-acetate (pH 7.9 at 25°C)  
50 mM potassium acetate  
10 mM magnesium acetate  
100 µg/ml BSA

### Unit definition

One unit is defined as the amount of enzyme required for complete digestion of 1 µg bacteriophage λ at 37°C for 1 hr in 50 µl reaction mixtures.

### Quality control

- Unit definition assay
- Overdigestion assay
- Endonuclease assay
- Extreme pure assay

### Dilution buffer:

FastGene® Diluent A

### Heat Inactivation

No.

### Methylation sensitivity

*dam* methylation: Not sensitive  
*dcm* methylation: Not sensitive  
CpG methylation: Not sensitive

### Prolonged incubation

A minimum amount of enzyme required to digest 1 µg substrate DNA for 16 hr; 0.5 U.

### Relative activity in FastGene® Buffers

FastGene® Buffer I: 75%  
FastGene® Buffer II: 100%  
FastGene® Buffer III: 75%  
FastGene® Buffer IV: 100%  
FastGene® FastCut Buffer: 100%

### Note

It is an isoschizomer of Hpa II. It is not affected by *dam*, *dcm*, or mammalian CpG methylation. However MspI can not digestion, when the external C in the sequence CCGG is methylated.

### Standard reaction condition

- Normal protocol

Component	Final Conc.	Volume
Substrate DNA	1 µg	X µl
10X FastGene® Buffer IV	1 X	5 µl
Msp I	20 unit	1 µl
Sterile water		up to 50 µl

→ Incubate at 37°C for 1 hr

- Fast protocol

Component	Final Conc.	Volume
Substrate DNA	1 µg	X µl
10X FastGene® FastCut Buffer	1 X	5 µl
Msp I	20 unit	1 µl
Sterile water		up to 50 µl

→ Incubate at 37°C for 15 min

※'Standard conditions' is only a general recommendation. The experimental conditions should be adjusted according to the purpose and sample.