

• **Software Key**

The software key comes with *Qsep* series instrument. Based on you order, you may get Basic or Advanced key.

Basic software key (blue string)



Advanced software key (green string)



Software key with different instrument type (*Qsep₁-Lite*, *Qsep₁*, *Qsep₁₀₀*, *Qsep₄₀₀*) tag has corresponding *Q-Analyzer* software.

	<i>Qsep₁-Lite</i>	<i>Qsep₁</i>	<i>Qsep₁₀₀</i>	<i>Qsep₄₀₀</i>
Tag				
Software				


Notification:

* *Q-Analyzer* and Drivers are for operating *Qsep* series instrument and must be installed before instrument connects to your computer.


*Each type of instrument can only be connected by its dedicated software.

1. Insert software key into your computer, enter the folder “Q-Analyzer” and double click the executable file to install the software.


* Please install the latest version of *Q-Analyzer* and the file name in this figure may different in your software key.

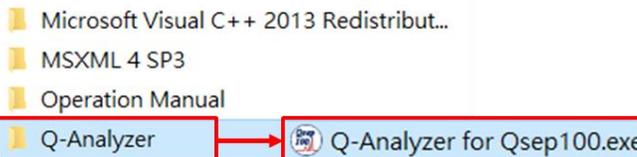


2. Select the folder to install *Q-Analyzer* (default folder is recommended)

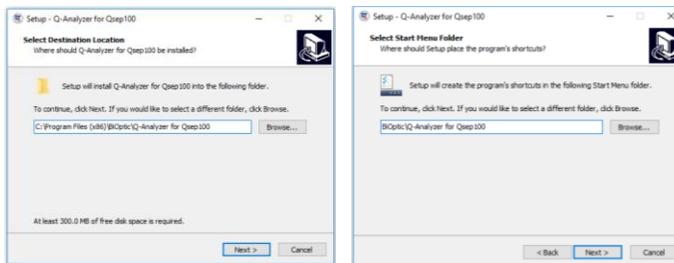

3. Select the components you want to install into the computer (Default settings are recommended to make sure *Qsep* instrument can operate smoothly).

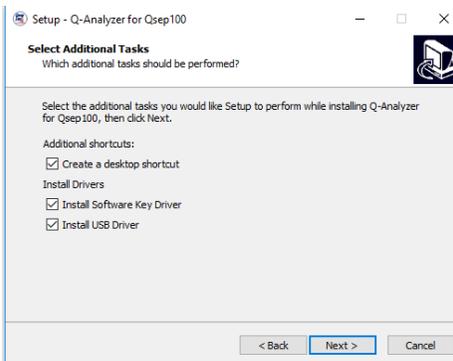


4. Confirm the settings and start to install *Q-Analyzer*.

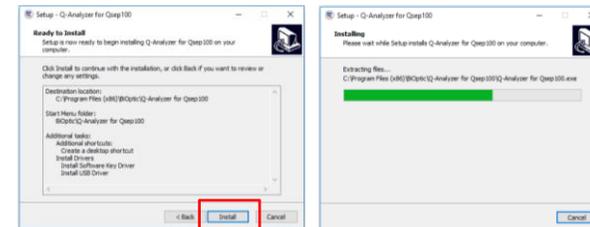

5. Start to install the drivers.



6. Please follow the instructions to complete the installation of Toolbox (Default settings are recommended).

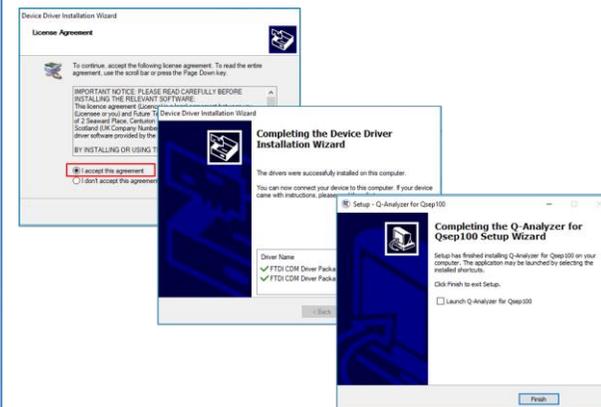


7A. Open *Q-Analyzer* without software key (Run as *Q-Analyzer* "Basic")

* Please make sure that "*Qsep* instrument" is connected to the computer.




7A-1. Double click "Q-Analyzer" icon on the desktop.

7A-2. Use *Q-Analyzer* "Basic" to operate *Qsep* instrument or analyze your data (For detail instructions please refer to "Quick start – Operation" or operation manual).

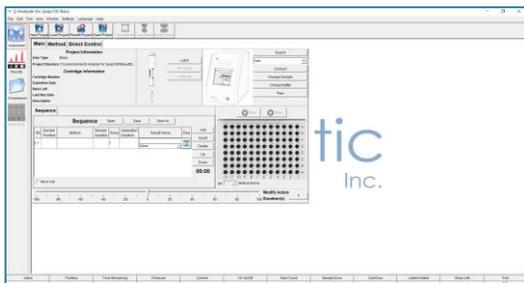

Notification:

This window will pop up when you try to start *Q-Analyzer* without software key or *Qsep* instrument connecting.

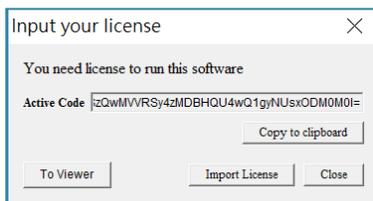

- You can import the license to enter *Q-Editor* or click "Close" to enter *Q-Viewer*.
- To activate "Advanced version" of *Q-Analyzer*, the software key need to be inserted into the computer.

7B. Open *Q-Analyzer* with software key (For *Q-Analyzer* "Basic" and "Advanced")


***Software key must be registered before use.**

Key Registration :



7B-1. Double click "Q-Analyzer" icon on the desktop.

7B-2. Using *Q-Analyzer* "WITH" software key, you need to fill in the blank in the registration window.

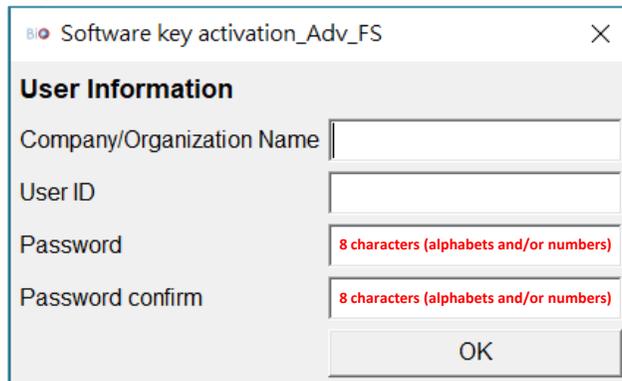


- Password must contain 8 characters, including alphabets and/or numbers
- The information will store at software key, not in the computer. You CANNOT change these information after pressing "OK".

If you have any question, please contact your local distributor or mail to "service@bioptic.com.tw" and ask for help.



7B-3. Double click *Q-Analyzer* icon on the desktop.

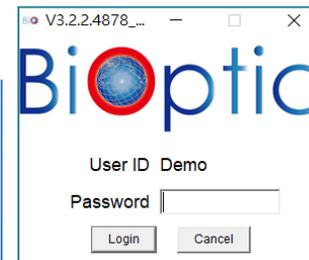


7B-4. Enter your password and start to use *Q-Analyzer*.

***WARNING: After Inputting wrong password 15 times, the software key will be locked.**

***If the software key is locked, please contact Bioptic**

7B-5. Use *Q-Analyzer* to operate *Qsep* instrument or analyze your data (For detail instructions please refer to "Quick start Operation" or operation manual).

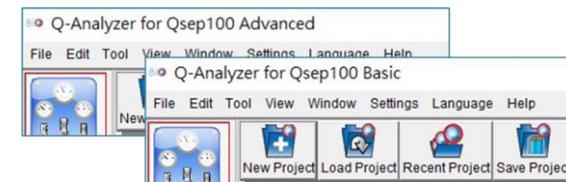

To use *Q-Analyzer* "Basic", you needs to connect at least one device below to your computer.



To use *Q-Analyzer* "Advanced", you **must** connect software key to your computer.



* Software key is required to activate *Q-Analyzer* Advanced.