

- *Q-Analyzer* installation includes *Q-Analyzer*, *Q-Editor* and *Q-Viewer*. The operation interface depends on your authority.
- *Q-Viewer* and *Q-Editor* are the software that allow user to read data (*Q-Viewer*) and to analyze data (*Q-Editor*) without software key or connecting *Qsep* instrument to the computer.
- *Q-Viewer* is a Freeware, but *Q-Editor* is NOT.

Basic software key (blue string)




Advanced software key (green string)



- **Software Key**
The software key comes with *Qsep* instrument.
* Based on you order, you may get Basic or Advanced key.

Q-Viewer


1. You can directly enter the *Q-Viewer* when the software key is not inserted and without connecting *Qsep* instrument.

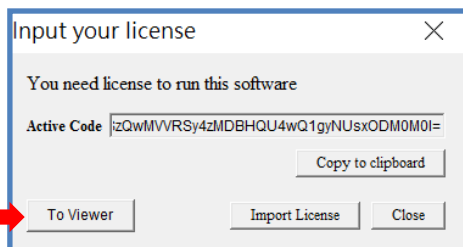


Q-Editor

The following steps are used for creating a license.

2. When you execute *Q-Analyzer* without software key and connecting *Qsep* instrument to the computer, "Input your license" window will pop out. Please click "Copy to clipboard".

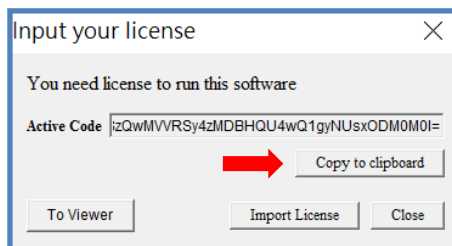


3. Insert the software key and click "Toolbox-License".

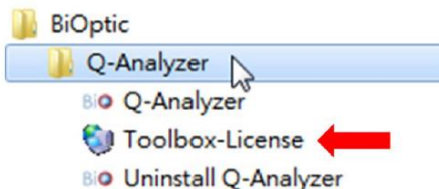


*"Toolbox" can be found in *Q-Analyzer* folder.
Default path: C:\Program Files (x86)\BiOptic\Q-Analyzer for Qsep


4. Click "Next".

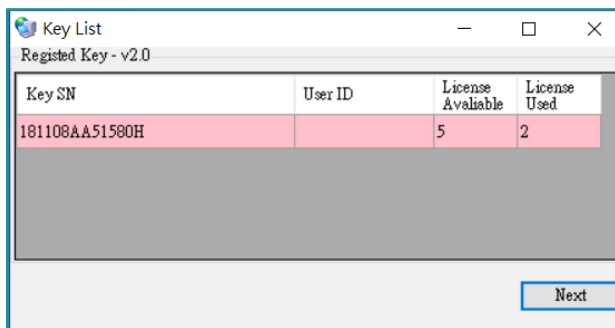


5. Paste the Activation Code (A). Fill in the "Remark" (B) and then click "Get New License" (C).

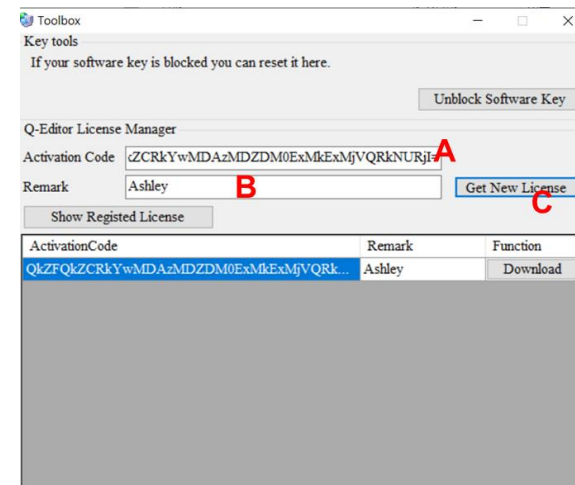


6. Enter the password of software key. If you forget the password, you can follow the unblock process. (Next page)

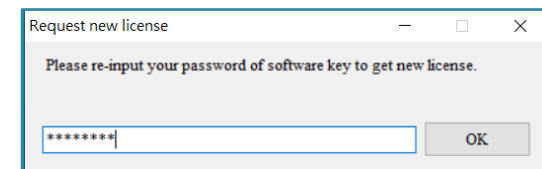


7. Once entering the code, the new license appears, please click "Download" to get the license.

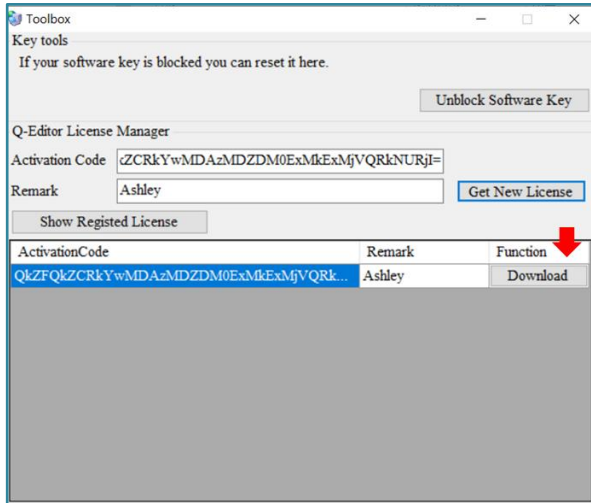

8. Back to the "Input your license" window, click "Import License", and then choose the license file you just downloaded.

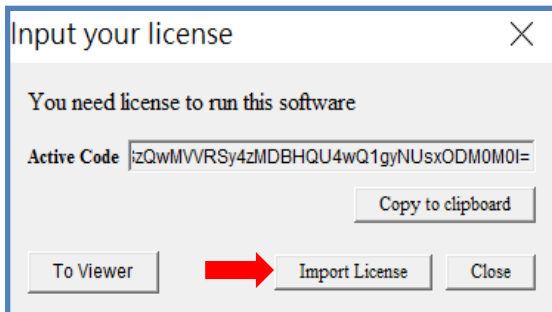


9. *Q-Editor* is ready to use.

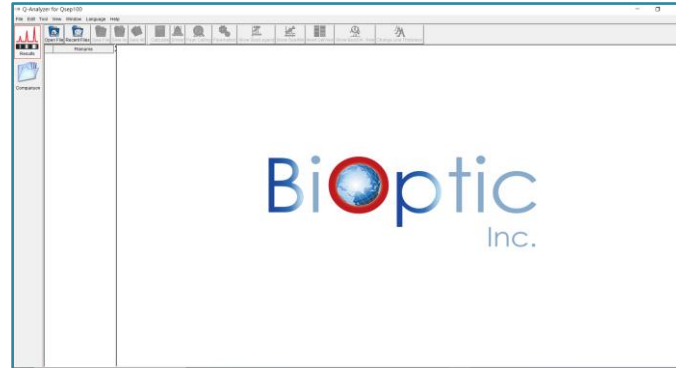


If you forget the password of software key, please follow the unblock process below.

Unblock process

- A. Click "Next".

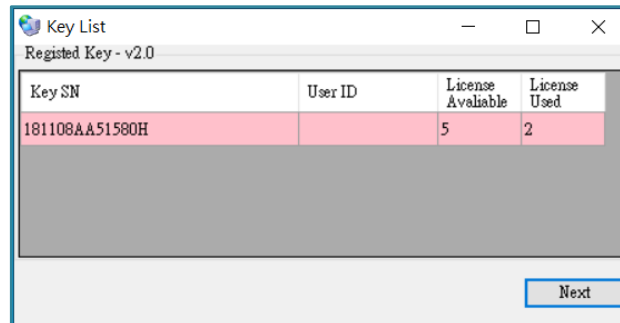

- B. Click "Unblock Software Key".

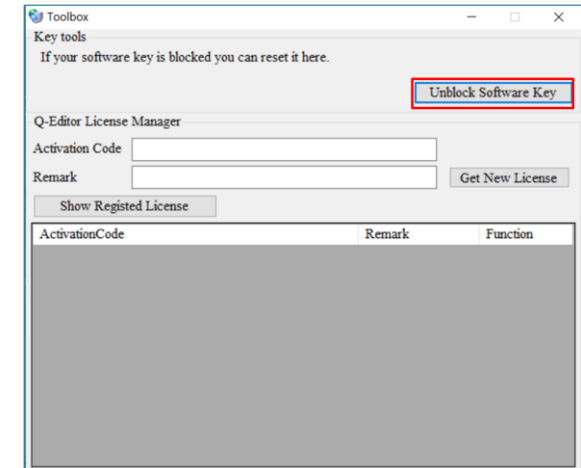

- C. The password will reset to "00000000".



Notification:

* Each software key provides **5 licenses**.
If you need more licenses of *Q-Editor*, please contact your local distributor.