


Update Procedure FAS-X 2.1.0

Important: It is recommended to perform a backup of the image files prior to performing an update.

- You need:
 - Empty USB drive in FAT32 format. Only use the USB-stick which is supplied with the FAS-X instrument for the update procedure.
 - Update File: **Update-FAS-X-2.1.0.zip**
- Unzip the **Update-FAS-X-2.1.0.zip** and copy the files onto the empty USB flash drive
- The files on the USB drive should look like this (Please note that the unzipped files contain another zip file: “Update.zip”. This should be left as it is):

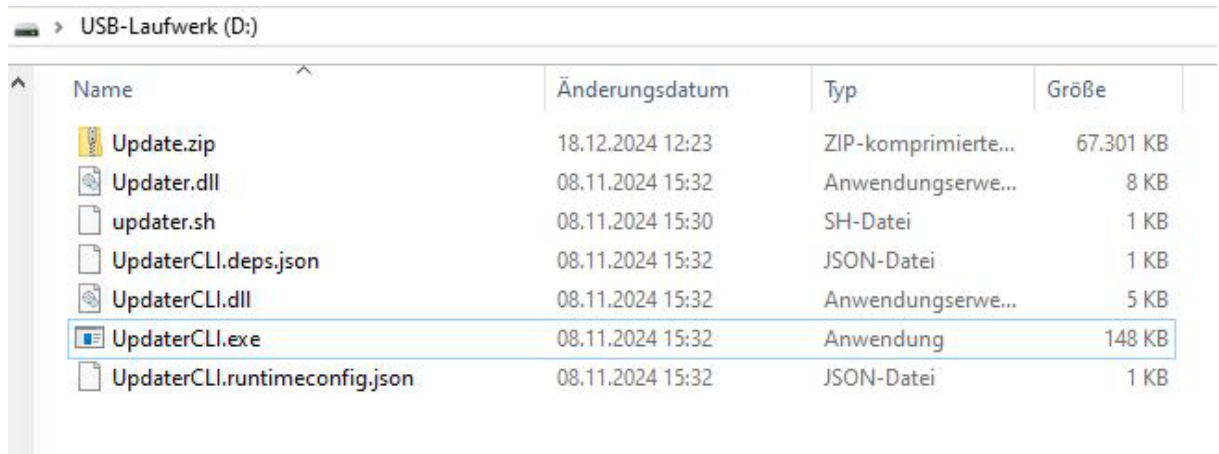

Name	Änderungsdatum	Typ	Größe
Update.zip	18.12.2024 12:23	ZIP-komprimierte...	67.301 KB
Updater.dll	08.11.2024 15:32	Anwendungserwe...	8 KB
updater.sh	08.11.2024 15:30	SH-Datei	1 KB
UpdaterCLI.deps.json	08.11.2024 15:32	JSON-Datei	1 KB
UpdaterCLI.dll	08.11.2024 15:32	Anwendungserwe...	5 KB
UpdaterCLI.exe	08.11.2024 15:32	Anwendung	148 KB
UpdaterCLI.runtimeconfig.json	08.11.2024 15:32	JSON-Datei	1 KB

- Start the FAS-X and Login as Admin
- Insert the USB Stick with the update files
- Go to **Preferences – Support – Admin – Managing**
- Press **Update Software**
- The FAS-X will update the software and reboot. This should take around 10-20 seconds.
- The new Software is now installed!
- Under **Preferences – Support** the **Version 2.1.0** should be displayed