


# FAS-X Image capture quick guide

## 3 taps for perfect gel image with Auto mode


- Connect the FAS-X to a power socket.
- Start the FAS-X software.
- Log in to your account. Once logged in, the software will open in Auto mode.  
*Green indicates the active mode.*




- Open the transilluminator drawer.
- Place your gel centrally.
- Close the transilluminator.
- Turn on the Blue/Green LED light.

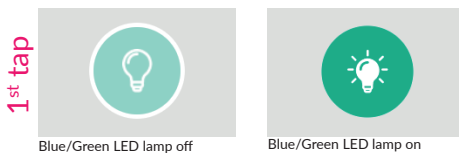

- Adjust the brightness.  
*5 different levels to adjust the brightness according to the signal intensity of the sample.*

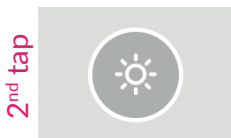

- Capture the image.

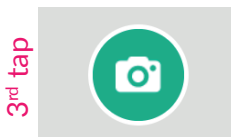

It's as simple as ABC!



## Get your personal demo of the FAS-X!

Get in touch with us and you will receive a complete product demonstration, or a demonstration adjusted to your specific needs!

[www.nippongenetics.eu/fas-x](http://www.nippongenetics.eu/fas-x)



## Access the complete FAS-X manual



<https://www.nippongenetics.eu/app/uploads/manual-gp-fas-x.pdf>

## Subscribe to the NIPPON Genetics EUROPE Newsletter!

Don't miss any news and stay informed about the latest product developments and current promotions.



[www.nippongenetics.eu/en/newsletter-registration/](http://www.nippongenetics.eu/en/newsletter-registration/)